

Title:	Shovel moves too far	ID:	
		<i>0191</i>	
Date in:	Response:	Model:	Author:
<i>2004-06-18</i>	<i>2004-06-18</i>	-	<i>CMa</i>

Q:

Quick question. Sometimes the shovel on our incubator serves the plate out a little too far onto the tray, so it's not sitting flat in its slot. Is there a way to adjust this, i.e. a data memory adjustment?

A:

The shovel is a hard-stop movement. It doesn't matter if the shovel moves out too far. In fact I prefer having the shovel come out too far since if it would move out not far enough it can lose the plate.