

| | | | |
|-----------------|---|---------------|----------------|
| Title: | StoreX Beeps after Sending the 1801 Commands | ID: | |
| | | 0212 | |
| Date in: | Response: | Model: | Author: |
| 2004-10-12 | 2004-10-12 | - | CMa |

Q:

In our driver we send a ST 1900 (Reset) and after we initialize (ST 1801) the StoreX will still beep. Any hints on this?

A:

The serial communication is interrupt driven. The microcontroller needs some time to update its internal flags – like the 1900-flag. In addition the 1900 command takes some time for execution. After you send a ST 1900 please add a delay (~0.1 sec.) then check the Ready (RD 1915). You may send the ST 1801 when the ST 1915 returns '1'.