

Title:	Carrousel Safety Switch through Door Interlock			ID:
				0027
Date in:	Response:	Model:	Author:	
2004-09-24	2004-09-24	Plate Hotel	CMa	

Q:

We are currently integrating an open air carousel into a work cell with guarding. It has been brought up that it would be useful to have the carousel by pass switch activate when the guarding door is opened. I would like to implement this using a interlock or limit switch. We want to ensure that we do not void warranty and we have your blessing when doing this modification. If this can be done I would require documentation on the carousel bypass switch to ensure the correct device is chosen.

A:

It's a standard NO (normally open) switch that switches 24V to Ground and a couple of Milliamps. An external Switch would have to connected parallel to this internal switch.