

Title:	Error 13 Hex			ID:
				0270
Date in:	Response:	Model:	Author:	
2007-05-02	2007-05-02	STX500	CMa	

Q:

Eric at AZ is having some problems with one of his older STX500's. Serial: 4003, FW ver: 6010. He has encountered the following error codes, when performing plate imports and exports:

6675 = 1A13h

6675 = 1A13h

Do you have any idea what could be the problem?

A:

Looks like the shovel hits something. Usually mechanical adjustment (shovel height, turn angle ..)