

|            |   |         |         |      |
|------------|---|---------|---------|------|
| Title:     | Deep Freezer Running at Plus Temperatures |         |         | ID:  |
|            |   |         |         | 0313 |
| Date in:   | Response:                                 | Model:  | Author: |      |
| 2008-10-21 | 2008-10-21                                | STX1000 | CMa     |      |

**Q:**

This STX1000 is currently running at -20C, we need this at +18C, I know you mentioned before we need to adjust the expansion valve, please send procedure ASAP so we can get our refrigeration guy to adjust this.

**A:**

No need for adjustment but check if there is no back-freezing at the compressor suction