

Title:	Lower Lift Position	ID:	
		0349	
Date in:	Response:	Model:	Author:
2010-02-19	2010-02-19	STX40 Shaker	CMa

Q:

We are trying to put 65mm plates into STX40 shaker Ser.Nr. 1524, it seems as we cannot go lower than $dm22=1$, I told him maybe move the zero reference lower, please read his response below.

A:

I assume this is a cabriolet type. The shovel motor height is over 60mm and has to fit under the open Gate. The lift should be able to handle this plate height.

However, if you'd to lower the lift height you may cut the black plastic block the lift initialized on. Beware of the spring force and possible interference if the lower end of the handler with the base plate.

