

Title:	<b>Database Backup</b>			ID:
				144
01.04.16	Response:	Model:	Author:	
		-	<i>CMa</i>	

## Database Backup

Q : Can you let me know what data needs to be backed up please. Is it a database, and if so does the database services need to be stopped before the backup. Or is it the contents of a folder that has normal files, and does anything have to be stopped before the data is backed up?

I am thinking of using a remote backup tool for each of the SAB PCs that we use for a lot of our servers, which would also benefit if a restore needs to be done, then it can be done directly to the PC.

A: The Database file is the only one that needs regular backup. The other files such as the software and log files may be backed up occasionally.

The database file can be copied during normal operation, the service does has to be stopped.

As I wrote in the response to Alex earlier this week there are two backup possibilities from our side:

- Automated task with defined frequency. We would need the network path and access rights to save files to this folder for backing up

- Place our database on a remote server which than can be backed up from your side. In this case network accessibility for our system to the database has to be ensured for proper functionality of the system.

All relevant data from our System can be found in the Drive D: on the PC. The database is in the Folder DB (Usually D:\Data\Liconic\DB).