
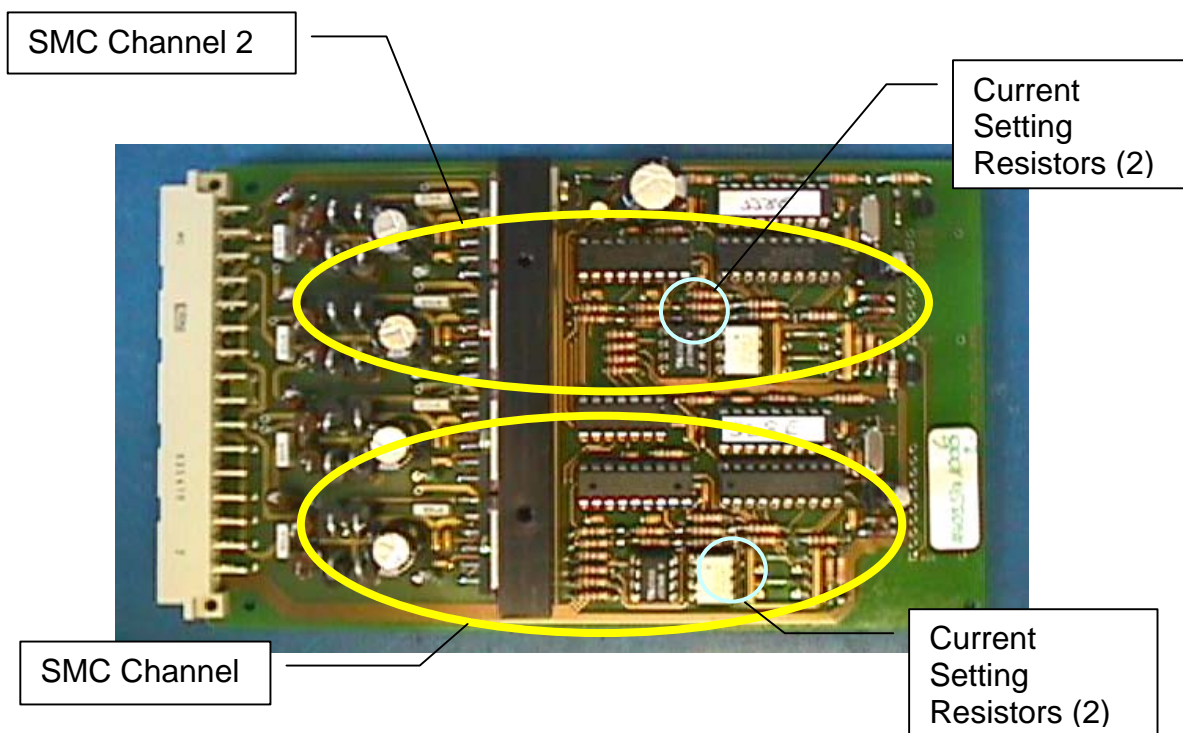


Title:	<b>SMC Motor Current Settings</b>			ID:
				0201
Date in:	Response:	Model:	Author:	
2004-08-20	2004-08-20	-	CMA	

## SMC Current Setting

The Motor Controller controls two axis (Channel 1 and Channel 2). The motor current is set for each channel by two resistors. Some cards have one axis (Channel 1) mounted only.

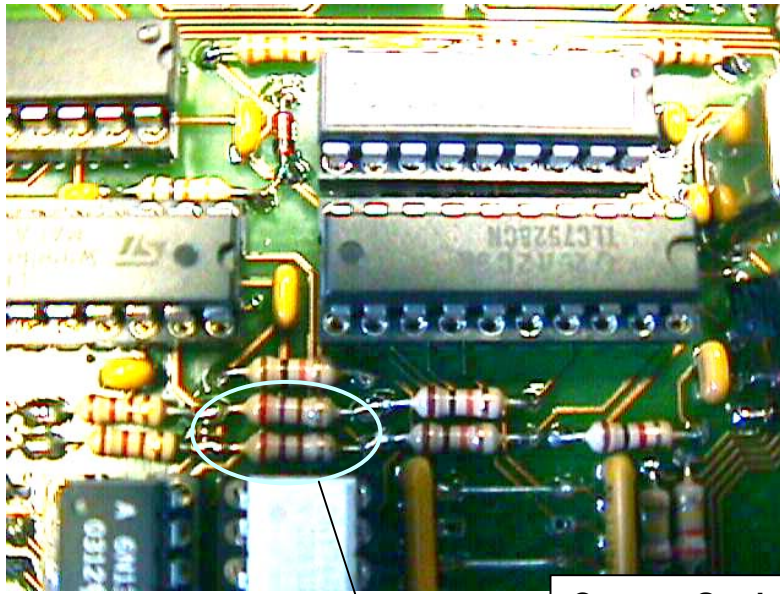

**Current Setting Resistors**

Low: 270R    red, purple, brown  
Med: 390R    orange, white, brown  
High: 680R    blue, grey, brown

To change the motor current remove the existing resistors and replace them with the needed value. Use Carbon or MOX resistors.