

<b>Title:</b>	<b>Overheating under the desk using STR</b>	<b>ID:</b>	
		0165	
<b>Date in:</b>	<b>Response:</b>	<b>Model:</b>	<b>Author:</b>
2006-02-15	2006-02-15	STR44	CMa

## Q:

We're having problems with an STR 44 HC overheating the EVO deck that it's underneath (see message thread below). Is there anything that can be done about it? this customer is going to purchase another 2 systems also, so as well as providing a retroactive solution we also need to take it into account when the next 2 are ordered. Please let me know if there's a solution for this problem.

## A:

I would add a big ventilator to get the heat out. I'm not familiar with the lower enclosure of the EVO. Make sure that you generate an airflow that pulls the generated heat out into the ambient.