

|            |                   |        |         |      |
|------------|-------------------|--------|---------|------|
| Title:     | STX Manual Access |        |         | ID:  |
|            |                   |        |         | 0225 |
| Date in:   | Response:         | Model: | Author: |      |
| 2004-04-05 | 2004-04-05        | STX200 | CMa     |      |

**Q:**

On the STX200 how do I manually turn the carousel to retrieve plates manually?

**A:**

Open the Front door(s) and rotate it manually. The motors are released when opening the glass door. You may also position the carousel by pressing the multi-function knob on the front panel. This will cause the carousel to move one position. If you press it for more than 1 second the carousel will rotate 1/3<sup>rd</sup> of a revolution.