

Title:	Carrousel does not stop rotating	ID:	
		0023	
Date in:	Response:	Model:	Author:
2003-09-19	2003-09-19	-	CMa

Q:

Unbeknownst to me, we had a 9 cassette StoreX unit in our loading dock. I just connected it and sent the initialization command. The carousel homes quickly, then slows down when it approaches cassette 1, but never stops!
Is there something wrong with this unit? Maybe the home sensor?

A:

No, you have to position the carrousel (e. g. "WR DM0 1") to stop it. The slow rotation is a leftover form the incubators that rotate by default.

If you want to check the sensor: It is an optical sensor that sits under the turntable. The PLC inputs are 000 and 001. You can read the sensor by "RD 0" or "RD 1". The sensor is on except if it "looks" through one of the holes of the turntable.