

Title:	Transferheight STR240	ID:	
		0306	
Date in:	Response:	Model:	Author:
2007-08-16	2007-08-16	STR240	CMa

Q:

You show the maximum reach of 1000, can this be changed using software?

A:

The maximum height we can reach is 1050mm (see also dimensions of the STR-Units). The height is variable by changing two data memories before initializing the unit.

- 1) DM22: This value defines the height of the Transfer station. Typical the value is 14500 (STR240). The absolute maximum for this value is 19000 (STR 240).
- 2) DM40: This is the value for the overdrive, which allows the Handling to reach over the top of the unit. Typical this value is 31000 (STR240). The absolute maximum for this value is 35500 (STR240).