

Title:	Climate Measurement Probe Tube on STX500 STX1000	ID:	
		0035	
Date in:	Response:	Model:	Author:
2006-03-04	2006-03-04	STX500 / STX1000	CMa

Q:

There is no hole in the glass door of my STX500 like there is on the STX40 or STX220. How can I take external measurement of the climate inside the incubation chamber?

A:

You may feed a tube between glass door and the front edge of the incubation chamber. Use a relatively stiff tube. We use PUR tube 2.5x5mm diameter.

