

Title:	DM24 DM22 Transfer Height does not work	ID:	
		0090	
Date in:	Response:	Model:	Author:
2005-10-10	2005-10-10	-	CMa

Q:

The value used for DM24 doesn't work.
I doesn't matter what we set the value to -- the shovel goes to the same height every time.
I've tried a whole range of values, only to arrive at the same result.

A:

Note that after a change of DM22 and/or DM24 you must initialize (ST 1801).
If the system is initialized already you must send an ST 1900 (Reset) and then an ST 1801
in order to make the system accept the new values.